

## Micropipette protocol card

For accurate measurements and to prevent damage to the micropipets, follow these important guidelines:

- Set the volume only within the range of your micropipet.
- Have the proper size disposable tip in place on your micropipet before immersion into any solution.
- Always keep the micropipet in a vertical position when there is liquid in the tip. In a horizontal position, fluid can leak back into the piston.
- Use your thumb to control the speed at which the plunger rises after taking up or ejecting liquid. Letting the plunger snap back damages the piston and the volume dispensed may be inaccurate.